V CARE CLEANCHECK

ENGLISH

LIQUID MEDIA FOR CLEANING CONTROL

INDICATIONS FOR USE

V CARE CLEANCHECK is intended for the control of cleaning procedures and for the training of operators.

PRODUCT DESCRIPTION

V CARE CLEANCHECK is a ready-to-use medium, which is fluorescent under UV-light (Black-Light).

COMPOSITION

- Ethanol 90%
- Non-Toxic fluorescent compound

PRODUCT SPECIFICATIONS

- Spectrophotometer test: passed
- Product appearance: passed
- A certificate of analysis and MSDS are available upon request.

WARNING

- Do not use after expiry date.
- Do not use the product if any signs of (microbial) contamination like decolourization, turbidity or a precipitate/cloudiness is observed.
- Do not use if the container is damaged upon delivery.

PRECAUTIONS

- Only for intended use.
- Always wear protective clothing
- The product may only be disposed of in accordance with the applicable regulations for waste products.
- Do not reuse solution that was already applied on surfaces

INSTRUCTIONS FOR USE

V CARE CLEANCHECK should be used only for its intended purpose by qualified and trained persons.

Control of cleaning procedures

- Open the vial of V CARE CLEANCHECK and dip the supplied swap in the media.
- Spread on the surface, which will be controlled for cleaning, before the cleaning procedure takes place.
- Wait at least 1 min fill the media, which was spread on the surface, is completely dry and not visible anymore.
- Conduct the regular cleaning procedure on the surface. *
- Use a UV-Light (black-light) source (we recommend a wavelength of 395 nm) to evaluate the surface. The media is fluorescent under UV-light. If no fluorescence is detected on the surface, the cleaning was conducted properly.

* For training purposes, the cleaning procedure can be conducted improperly to demonstrate the importance of following the procedure correctly.

STORAGE INSTRUCTIONS AND STABILITY

- Store between 2°C to 37°C.
- Keep away from sunlight.

FEEDBACK AND OBLIGATION TO INFORM

Thank you for purchasing our product. If you have any suggestions for us to improve our products, please send them to qc@vitromed.com.

Report any serious incidents that have occurred with the product immediately by email to qc@vitromed.com.

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VITROMED GmbH